IVIS Living Image download, installation, and License Seat Maintenance

Users can download Living Image 4.7.4 for FREE

For a permanent license of Living Image, you will need a 20-digit license key to activate this product - Contact SAIRF@uwcarbone.wisc.edu for the license key, and include your PI initials (IVIS login) to receive Research Drive network access to your data.

To download, use the link below. You must be logged in as the administrator to download this software (contact your IT Dept if you are not an admin). Choose the appropriate installer corresponding to your operating system. After the file is downloaded, right click on the file and select "Run as administrator." Choose the default settings and complete the installation.

https://www.perkinelmer.com/lab-products-and-services/resources/in-vivo-imaging-software-downloads.html#LivingImage

<u>Living Image 4.7.4 Windows 64-bit installer</u> - when installing, <u>do not</u> check the boxes for Acquisition

Living Image® Software (IVIS Imaging Systems)		
Living Image 4.8.0 (for use on all IVIS systems except IVIS Spectrum)		
Setup Files	User Guides	Release Notes
Windows x64 (64-bit)	Installation & Licensing Guide (PDF)*	4.8.0 Release Notes (PDF)
MacOS	ReadMe.txt	
Living Image 4.7.4 太		
Setup Files	User Guides	Release Notes
Windows x64 (64-bit)	Installation & Licensing Guide (PDF) ReadMe.txt	4.7.4 Release Notes (PDF)
Living Image 4.7.3		
Setup Files	User Guides	Release Notes
Windows x86 (32-bit)	Installation & Licensing Guide (PDF)	4.7.0 Release Notes (PDF)
MacOS	ReadMe.txt	4.7.1 Release Notes (PDF)
		4.7.2 Release Notes (PDF)
		4.7.3 Release Notes (PDF)
*LI 4.8.0 user guide coming soon. Please refer to LI 4.7.4 installation & licensing guide.		